Resident Case Distribution Beaker AP Guide:

Assigning the cases to different residents/fellows:

1. Click on your **Outstanding List** tab or find the link to the **Outstanding List** from the Beaker AP Lab Dashboard.



2. Filter the list so it only shows the Case Status of Gross Done:

	Case Status	Docidont 🔺	E Test Details
SYN	Accessioned	Ascending	- M rest Details
ES TU	Accessioned	✓ Descending	
ON P	Accessioned	Custom	SP18-00
EDLE	Accessioned	Filter by 🔹 🕨	(AII)
ES TU	Gross Done	Clear all filters	📕 (Blanks)
ON P	Accessioned		(NonBlanks)
ON P	Accessioned		Accessioned 2
	Gross Done		Amendment
ES TU	Accessioned		Gross Done er
Y	Accessioned	Todd P Willian	Path Rev Com
INS A	Gross Done	Todd P Willian	Res Reviewed
ECTO	Amendment	Homunculus E	

- a. Right click on the gray Case Status column header.
- b. Click on Filter By.
- c. Choose **Gross Done**. This will hide all the other cases so only the cases with the Gross Done status will be showing.
- 3. Once you have limited the Outstanding List to only show the Case Status of "**Gross Done**" you can select multiple cases at a time.
 - a. To select several cases that next to each other, click on the first case and hold down the SHIFT button on your keyboard. Click on the last one. This should select all the cases in between those two cases.

SP17-00840	Beaker, John	A - LIVER BIOPSY	Gross Done		
SP17-00845	Beaker, Mimi	A - LIVER BIOPSY, B - E	Gross Done		
SP17-00847	Beaker, Bubba	A - LIVER BIOPSY, B - LI	Gross Done		Susan Woods
SP17-00864	Beaker, Mama Walo	A - KIDNEY ALLOGRAFT	Gross Done		Rita M Braziel, MD
SP17-00867	Kidney, Augtest A	B - APPENDIX INCIDENT	Gross Done		Rita M Braziel, MD
SP17-00876	Beaker, Corelle	A - LIVER BIOPSY, B - B	Gross Done		Rita M Braziel, MD
SP17-00885	Beaker, Alien	A - APPENDIX INCIDENT	Gross Done		
SP18-00058	Beaker, Jenny	A - VAGINAL BIOPSY	Gross Done		Amanda K Vansand
SP18-00054	Transplant Test, Kri	A - BRAIN BIOPSY	Gross Done		Marjorie R Grafe, M
ID	Name	Protocols	Case Status	Resident	Pathologist

b. To select different individual cases, hold down the CTRL key and single click on each case you want to select.

ID 🛕	Name	Protocols	Case Status •	Resident	Pathologist
SP17-00834	Beaker, Alicia	A - LIVER ALLOGRAFT B	Gross Done	Homunculus Beaker	Herbert West Beaker
SP17-00835	Beaker, Jenny	A - LIVER BIOPSY, B - LI	Gross Done	Homunculus Beaker	Herbert West Beaker
SP17-00836	Beaker, John	A - LIVER BIOPSY	Gross Done		
SP17-00837	Beaker, John	A - LUNG WEDGE BIOPSY	Gross Done	Todd P Williams, MD	
SP17-00838	Caspian, Sea	C - BREAST MARGINS A	Gross Done	Todd P Williams, MD	Daniel T Fitzpatrick
SP17-00839	Beaker, John	A - LIVER BIOPSY	Gross Done		Donald C Houghton,
SP17-00840	Beaker, John	A - LIVER BIOPSY	Gross Done		
SP17-00845	Beaker, Mimi	A - LIVER BIOPSY, B - E	Gross Done		
SP17-00847	Beaker, Bubba	A - LIVER BIOPSY, B - LI	Gross Done		Susan Woods
SP17-00856	Beaker, Harold	A - BREAST BIOPSY	Gross Done	Todd P Williams, MD	

4. Once you have selected the cases, click on the triangle next to the **L** Assign button and pick Assign.

1 Assign - D
🛉 <u>T</u> ake
Put Back
💄 Assign

5. Choose the resident/fellow's name and press **Assign** button.

Assign Case - SP17-00847				
<u>U</u> ser:	WILLIAMS MD, TODI	ļ ļ		
<u>R</u> ole:	Resident Pathologist	Resident Pathologist 🔎		
	Assign <u>C</u> a	ancel		

6. Continue assigning cases until all the cases have been assigned.

© 2017 Epic Systems Corporation. All rights reserved. PROPRIETARY INFORMATION - This item and its contents may not be accessed, used, modified, reproduced, performed, displayed, distributed or disclosed unless and only to the extent expressly authorized by an agreement with Epic. This item is a Commercial Item, as that term is defined at 48 C.F.R. Sec. 2.101. It contains trade secrets and commercial information that are confidential, privileged and exempt from disclosure under the Freedom of Information Act and prohibited from disclosure under the Trade Secrets Act. After Visit Summary, Analyst, App Orchard, ASAP, Beaker, BedTime, Bones, Break-the-Glass, Caboodle, Cadence, Canto, Care Everywhere, Charge Router, Chronicles, Clarity, Cogito ergo sum, Cohort, Colleague, Community Connect, Cupid, Epic, EpicCare, EpicCare Link, Epicenter, Epic Earth, EpicLink, EpicWeb, Good Better Best, Grand Central, Haiku, Happy Together, Healthy People, Healthy Planet, Hyperspace, Identity, IntraConnect, Kaleidoscope, Limerick, Lucy, MyChart, OpTime, OutReach, Patients Like Mine, Phoenix, Powered by Epic, Prelude, Radar, RedAlert, Resolute, Revenue Guardian, Rover, SmartForms, Sonnet, Stork, Tapestry, Trove, Welcome, Willow, Wisdom, and With the Patient at the Heart are registered trademarks, trademarks or service marks of Epic Systems Corporation in the United States of America and/or other countries. Other company, product and service names referenced herein may be trademarks or service marks of their respective owners. U.S. and international patents issued and pending.

This guide is based on Epic 2017.